

Please do not remove the pre-existing thermometer and probes until the new thermometer is confirmed to be functioning properly.

TraceableLIVE® BASIC SET UP INSTRUCTIONS


DEVICE SET UP

1. PROBE CONNECTION: Connect “P1” probe into “P1” slot and “P2” into “P2” slot on the top of device (serial numbers on the probes must match the serial number on the unit)



2. BATTERY ACTIVATION: Remove the 2 plastic slips from the battery compartment to activate the batteries.



3. VIEW THE TEMPERATURE FOR THE APPROPRIATE PROBE CHANNEL: Press Channel Select  to display the desired channel, or dual channels.

4. SELECTING THE UNIT OF TEMPERATURE MEASURE: Press **C/F** select the desired unit of temperature measure (°C or °F).



1 WiFi Setup



Press/HOLD WiFi button "AP" will appear on display.

Press WiFi button again, "AP URI" will appear on display.

When "AP rEAdY" appears, proceed to next step.

Back of unit

EXAMPLE

6500CC
S/N 100080013
KEY 5555123456
.....
traceable.com 281-482-1714
By: 100 Due: Nov, 19 2024

EXAMPLE

2 Use Internet enabled device to connect to WiFi



Select Network:
CC6500-0013
see example network

Example
Network: CC6500-0013

EXAMPLE

6500CC
S/N 100080013
KEY 5555123456
.....
traceable.com 281-482-1714
By: 100 Due: Nov, 19 2024

EXAMPLE

3 Open Browser, type 192.168.1.1 in address bar

Select local network and enter password.
Security type WPA2 recommended or contact your network admin.



4 Login to your Account traceablelive.com



CONNECT DEVICE TO WIFI AND ACCOUNT LOGIN

Device will restart.

DISPOSAL OF PRE-EXISTING THERMOMETER: After confirming the new thermometer is functioning properly, the pre-existing thermometer and probes can be disposed:

- Put old thermometer in zip seal bag used for hazardous waste
- Place the bag in the "Rx" Hazardous Waste tote

Need additional Assistance?
Cole-Parmer Houston Customer Service
Phone: 281-482-1714 from 8:30 am to 5:00 pm CST
Email: support@traceable.com